


MODEL M45 PUMP

Installation and Maintenance Instructions

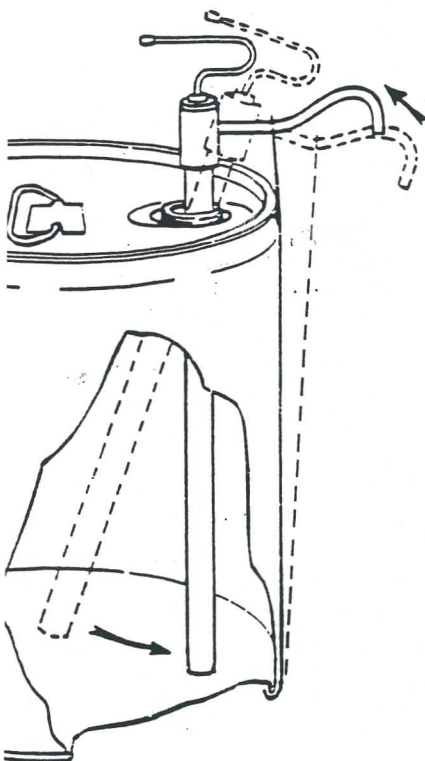

1. Slip the band over spigot of pump, making certain that the bottom bend on band faces pump.
2. Insert pump barrel in bail opening at angle
3. When used with RIEKE® FLEXSPOUT® closure, screw black cap provided on pump cylinder to closure for secure fit.
4. Slip end of band under bottom chime of pail.
5. Holding spout of pump, straighten pump to upright position to tighten band.
6. Packing nut at shaft may be tightened when required to prevent leakage.
7. If pump does not prime quickly, take the following corrective measures:
 - a. Rinse with hot water to remove possible solidified deposits in check valve.
 - b. Disassemble pump, inspect plastic piston for damage, and check to be sure ball in foot valve is operating freely.
8. Plastic spout cap prevents evaporation and residual drippage. Replace cap after dispensing desired amount of material.
9. If pump is removed from pail, rinse thoroughly in hot water before again placing in use.
10. CAUTION: Metal pumps are not recommended for use with acid products.