

HOW TO ASSEMBLE THE DRYING RACK 213

After inserting the expansion dowels, assemble the drying rack as directed herebelow:

- 1) Fit the drain pipe connection (see picture **A**).
- 2) Insert the drying rack into the lower and upper dowels, then screw tight the fastening knob to the lower expansion dowel (see picture **B**).
- 3) Insert pegs to complete assembly (see picture **C**).
- 4) In case you need two drying racks side by side, assemble them (after performing step 1) and couple them up by using the accompanying fastening stopper (see picture **D**).

Note: Drill Template for fixing the Drying Rack to the wall is included in the Drying Rack Packaging.

