

Kartell LABWARE



Pipette & Burette Rinsing Set

www.kartell.com.au

Pipette & Burette Rinsing Set

- Designed for washing pipettes & burettes up to 600mm long.
- Range of sizes available to suit the application.
- Polyethylene rinsers (Art. 215 & Art. 217) fill via the spray nozzle situated on the top rim, eliminating back-syphoning as the water inlet is located above and away from any contaminated waters.
- All contaminated water syphons out of the rinser through the specially designed syphon pipe leading to the large bore outlet tube situated underneath the rinser.
- The rinser refills automatically during this process.
- Pipettes and burettes are contained in specially designed baskets during washing to minimise breakage.

The washing process is as follows:

- Place the pipettes in the basket (*Art. 219 or Art. 222*).
- Place the basket into the jar (*Art. 218*) full of washing solution. Leave to soak for 2-3 hours so the solution can clean the pipettes.
- Remove the basket from the jar (*Art. 218*) and place in the rinser (*Art. 215 or Art. 217*). Connect the water pipe to the spray nozzle, and the outlet tube to a drainage point.
- Fill the rinser with water, add detergent solution and shake several times to ensure a thorough washing.
- Turn the water tap on and allow rinsing cycle to complete.

Full instructions are included with each product in the range, containing more details depending on the application required.

Suggested sets consist of the following products:

Pipettes – 1 x 217, 2 x 219 and 1 x 218

Burettes – 1 x 215, 2 x 219, 2 x 1219 and 1 x 218

PIPETTE & BURETTE RINSING SET				
art.	Description	Diameter Ø (mm)	Basket height (mm)	Height (mm)
215	Automatic Burette Washer	170	-	990
217	Automatic Pipette Washer	170	-	734
218	Pipette Jar	162	-	650
219	Pipette Basket - Long Handle	145	300	648
221	Pipette Jar	162	-	503
222	Pipette Basket	145	300	497
223	Pipette Jar	125	-	250
1219	Extension Handle for Art. 219	-	-	290

For more information on these products, please see page 74 of the 2012/2013 KARTELL catalogue, available for download at:

www.kartell.com.au



Available From:



SIEPER Group
Wholesale Distributors

ABN: 58 000 296 453

101 - 109 Deakin St Silverwater NSW 2128
 PO Box 6724 Silverwater NSW 1811
 (02) 9352 0600 (02) 9748 0566
 sales@sieper.com.au www.sieper.com.au